



## NeutraGreen™

Liquid Aldehyde Neutralizer  
Instructions for Use

[statlab.com](http://statlab.com)

NeutraGreen™ neutralizes waste formalin to a non-hazardous material. The formaldehyde solution is irreversibly changed to a non-hazardous material and cannot change back to formaldehyde. Treated waste formalin is odorless. NeutraGreen™ neutralizes formaldehyde to 1% concentration in two hours and is completely neutralized after 4 hours. Environmental testing proves NeutraGreen™ will neutralize waste formalin to not only a non-hazardous but a non-toxic material as well.

### APPLICATIONS

- 10% Neutral Buffered Formalin
- 20% Neutral Buffered Formalin\*
- 10% Glutaraldehyde Solution\*

\*may require additional time for complete neutralization

### NeutraGreen™ PACKAGING

NeutraGreen™	1525-18	5 Gallon Cubitainer
NeutraGreen™	1525-25	2.5 Gallon Cubitainer
NeutraGreen™	1525-41	4 x 1 Gallon Bottles

### INSTRUCTIONS FOR USE

1. Pour waste formalin into a sturdy non-leaking container with enough headroom to add NeutraGreen™ at a ratio of 1 part NeutraGreen™ to 4 parts formalin.
2. Add NeutraGreen™ in the proper ratio.
3. Wait 2 hours and test for neutralization. If less than 1% aldehyde present product is ready to be discarded. All formalin vapors should be noticeably gone.
4. If desired, it is recommended to perform testing to ensure that product has been neutralized.
5. Once formalin is neutralized, material may be disposed of as non-hazardous and non-toxic material with zero vapor exposure. Dispose of waste material according to all local, state and federal rules and laws.

Technical Note: dispose down sanitary sewer if permitted.

### STORAGE/ SAFETY

Storage: Room Temperature

Refer to SDS for details-Always wear chemical resistant gloves and eye protection when handling chemicals

Follow all local, state and federal laws and regulations regarding formalin solutions.

### ORDERING INFORMATION

Product	Product Number	
NeutraGreen™	5 Gallon Cubitainer	1525-18
NeutraGreen™	2.5 Gallon Cubitainer	1525-25
NeutraGreen™	4x1 Gallon Bottles	1525-41

### FORMALIN NEUTRALIZATION TEST PRODUCTS AVAILABLE FROM STATLAB:

2183100 Formaldehyde Test Kit (NeutraGreen™)  
(Range 0.05-1% or 0.5-10%)

V1.0

### CONTACT INFORMATION

StatLab Medical Products, Inc.  
407 Interchange Street  
McKinney, Texas 75071  
Phone: 800-442-3573 Fax: 972-436-1369  
[www.statlab.com](http://www.statlab.com)

Should any of our materials fail to perform to our specifications, we will be pleased to provide replacements or give credit for the purchase price. The information given in this bulletin is to the best of our knowledge accurate, but no warranty is expressed or implied. It is the user's responsibility to determine the suitability for their own use of the products described herein, and since conditions of use are beyond our control, we disclaim all liability with respect to the use of any material supplied by us. Nothing contained herein shall be construed as a recommendation to use any product or to practice any violation of any law or government regulation.

2015-March 12