

OdorVaultTM

Lock Up Those Odors

(Rev. 12/6/18)




ODORVAULT™, THE PRODUCT AND ITS PURPOSE:

OdorVault™ is a revolutionary new product that is designed to absorb the worst odors! The OdorVault™ can be placed onto or near any surface, such as: trash cans, bio-hazard bins, fume hoods, cabinets, shelves, and refrigerators, etc. There are three options to choose from when purchasing the OdorVault™. The OdorVault™ Formalin, which is great for formalin odors, the OdorVault™ Xylene for xylene vapors, or the OdorVault™ Organic, which is perfect for all other foul odors.


ODORVAULT™, INSTRUCTIONS FOR USE:

The **OdorVault™** can be used in two ways: with a box or with an adhesive strip that can be attached to any surface. Make sure to clean the surface where the **OdorVault™** will be placed.

OPTION #1: ODORVAULT™ BOX:

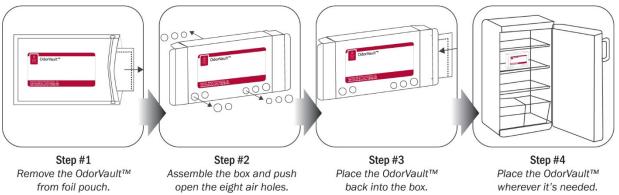

OPTION #2 ODORVAULT™ ADHESIVE STRIP:

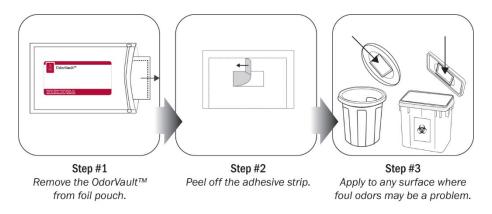

ODORVAULT™, PRECAUTIONS:

Store the **OdorVault**[™] at room temperature away from direct sunlight. Do not open the contents of the **OdorVault**[™] bag. **OdorVault**[™] can't be used in areas that are wet, damp, or exposed to high temperatures/high humidity. The minimum temperature to which the **OdorVault**[™] Adhesive Strip can attach to is 10°C. The adhesive strip strength may be weakened by excessive re-positioning or hand contact. Do not adhere to objects that may break easily.

Copyright ©2012-2019 American MasterTech Scientific. All rights reserved.

No part of these pages may be used for any purpose other than personal use. Therefore, reproduction, modification, storage in a retrieval system or retransmission, in any form or by any means, for reasons other than personal use, is strictly prohibited without prior written permission.

ODORVAULT™, DISPOSAL:

Peel away the **OdorVault™** slowly or remove from the box and place it back into the foil pouch that it came in. Discard it as industrial waste. If adhesive residue remains on the surface, wipe it off with an adhesive cleaner.

ODORVAULT™, **STABILITY**:

The **OdorVault™** will last up to three months.

CONTACT INFORMATION:

American MasterTech Scientific 1330 Thurman St. Lodi, CA 95240 USA Tel 800 860 4073 Fax 209 368 4136 www.americanmastertech.com StatLab 2090 Commerce Drive McKinney, TX 75069 USA Tel 800 442 3573 Fax 972 436 1369 www.statlab.com