



Treated Waste Aldehyde Test Kit Instructions

1. Draw sample of treated aldehyde waste with pipette and fill square mixing bottle provided in kit to the **10ml** mark.
2. Open one (1) Aldehyde Reagent Powder Packet and pour contents into the square mixing bottle.
3. Secure the mixing bottle opening with glove-thumb or cap and swirl to mix solution for 10 seconds.
4. Add two (2) drops of color indication solution (thymolphthalein) to the mixing bottle, swirl gently and note color.
5. If aldehyde is present the sample will turn BLUE.
6. If sample turns blue, secure the lid on the FRC-3™ collection container, agitate solution and allow solution to stand for additional 20 minutes, and then re-test.
7. If sample does not turn blue, then aldehyde waste solution is now ready for discharge into the sanitary sewer.