

MAINTENANCE

1. Always unplug Paraffin Dispenser before moving or cleaning. HAZARD – Having water or any other liquid near or in contact with power supply may cause electrical shock
2. Empty and clean Paraffin Dispenser regularly. To clean use a lint-free type of towel and sponge when necessary. Avoid using any type of instrument/tool/chemical that can scratch or damaged the Paraffin Dispenser.
3. The Paraffin Dispenser has been carefully tested and inspected.

WARRANTY

The manufacturer warrants this instrument to be free from defects in material and workmanship under normal use for one year from the date of purchase. It does not cover damage resulting from abuse or misuse, repairs or alterations performed by other than authorized repair technicians, or damage occurring in transit. If you have questions concerning this product or warranty, contact the dealer from whom it was purchased. For warranty service, Paraffin Dispenser should be well packed to avoid damage in transit, preferably in original box and packing. Include your complete return address and telephone number as well as a description of the difficulty, date and place of purchase, and ship to the address below. It will be repaired or replaced at no charge and returned. If misuse, alterations, accident or abnormal conditions of operation caused failure, an estimate for repairs will be provided for your approval prior to work being performed.

Premiere Service Department
7241 Gabe Court
Manassas, VA 20109
(703) 330-1413

Premiere®



XH-4002, 2.3 Gallon Paraffin Dispenser

USER'S MANUAL

Before using this product, please read this manual and learn about all the available features to obtain the best results from your equipment.

Warning:

To prevent fire or shock hazard, do not expose this unit to rain or moisture. This equipment should be used with AC 110V, 60Hz in the USA or Canada.

No user-serviceable parts inside. Refer servicing to qualified service personnel.

Premiere® trademark reg. No. 1713212 USA

APPLICATION

This Paraffin Dispenser is designed for melting paraffin and dispensing the paraffin. The sole purpose of this Paraffin Dispenser is for histology use along with research, instruction, and experiments in universities and technical secondary schools. The Paraffin Dispenser is not recommended for any other uses than the ones described below in this user manual


ASSEMBLY INSTRUCTIONS

1. Remove Paraffin Dispenser from Styrofoam packing and place it on a stable worktable. Remove all plastic bags and paper covering (these can be discarded).
2. Connect the power supply cord to the unit and to an 110V outlet and your **Premiere**® Paraffin Dispenser is ready for use.

OPERATION

1. Prior to use fill the Paraffin Dispenser with solid paraffin at least halfway full. It is also recommended not to fill the Paraffin Dispenser with too much paraffin so that it has the chance to overflow. **HAZARD** – Having liquids near or in contact with power supply may cause electrical shock.
2. After making sure that the Paraffin Dispenser is plugged in, turn on the machine with the switch on the front. The switch indicator light (#3) should light up to indicate the machine is working.
3. Before using check the temperature to make sure that your set temperature is above the melting point of your paraffin (as indicated on the paraffin box.) A thermometer can be inserted in the hole in the lid.
4. To adjust the temperature, press the set button (#4). Then use the up and down arrow buttons (#5) to set the temperature at the desired setting. This adjustment should be visible in the setting display panel (#2). The actual temperature of the wax basin can be seen in the actual temperature display panel (#1). Please allow ample time for the Paraffin Dispenser to warm up and start melting the paraffin.
5. Once the Paraffin Dispenser reaches your set temperature, the heat light indicator (#3) will go out.

6. This Paraffin Dispenser is also built with a timer range. You can set the dispenser to run for a certain amount of time. Once again if you push the set button (#4) the setting display panel should change to a timer screen (#2). Then use the up and down arrow buttons (#5) to set the timer at the desired setting. This adjustment should be visible in the setting display panel (#2). Once you reached the desired time, hit the set button (#4). Your timer will count down and buzz (meaning it turned off) when your time has elapsed.

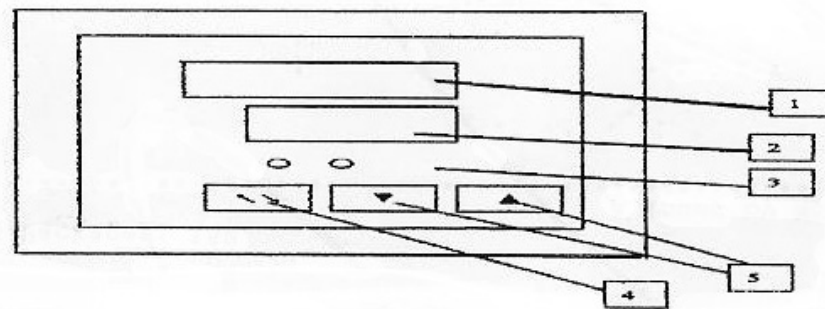

- 1) Temperature Display
- 2) Setting Display
- 3) Switch Light Indicator
- 4) Set Button
- 5) Up and Down arrow buttons

Cautions

- 1) Unplug the Paraffin Dispenser when not in use.
- 2) Do not use the faucet when the temperature is below the melting point of the paraffin.
- 3) Be very careful with the faucet, as it may be hot because it can absorb the heat from the Paraffin Dispenser when the dispenser is set at high temperatures.
- 4) Please contact your dealer with any operating problems. Only authorized repair technicians should perform repairs.